20190925
Ver. 0 (KR)

AccuPrep[®]

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8-11, Munpyeongseoro, Daedeok-gu, Daejeon, 34302, Republic of Korea
Tel: (Korea) 1588-9788 (International) +82-42-930-8777 Email: sales@bioneer.com

BIONEER
Innovation • Value • Discovery

AccuPrep® Kit 와 함께

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Selection Chart

Nucleic Acid	Sample	Product Type	Product	Yield	Experiment Time
Genomic DNA	Cultured animal cells Lymphocytes Mammalian tissue Whole human blood and bacterial cultures	Spin Column	AccuPrep® Genomic DNA Extraction Kit	Up to 20 µg	< 30 min
	Plant tissues	Spin Column	AccuPrep® Plant Genomic DNA Extraction Kit	Variable*	< 50 min
	Fresh or frozen stool Samples with large amounts of PCR inhibitors	Spin Column	AccuPrep® Stool DNA Extraction Kit	2 - 5 µg	< 50 min
	Cultured animal cells Lymphocytes Mammalian tissue Whole human blood and bacterial cultures	96-Well Plate	AccuPrep® Genomic DNA Extraction Kit for 96 well vacuum Block	Variable*	< 40 min
Plasmid DNA	Bacterial cultures	Spin Column	AccuPrep® Nano-Plus Plasmid Mini Extraction Kit	Up to 30 µg	< 10 min
		Spin Column	AccuPrep® Nano-Plus Plasmid Midi Extraction Kit	Up to 75 - 100 µg	< 40 min
		Pump			< 30 min
		Spin Column	AccuPrep® Nano-Plus Plasmid Maxi Extraction Kit	Up to 300 - 500 µg	< 60 min
		Pump			< 45 min
		Spin Column	AccuPrep® Plasmid Mini Extraction Kit	Up to 20 µg	< 20 min
96-Well Plate	AccuPrep® Plasmid Mini Extraction Kit for 96 well vacuum Block	Up to 10 µg	< 40 min		
Fragment DNA	PCR products DNA fragments in agarose gel slice Labeled, modified or digested DNA	Spin Column	AccuPrep® PCR/Gel DNA Purification Kit	70 - 90% recovery	< 30 min
	PCR products Labeled, modified or digested DNA	96-Well Plate	AccuPrep® PCR Purification Kit for 96 well vacuum Block	70 - 90% recovery	< 25 min
RNA	Cultured animal cells Mammalian tissue Plant tissue	Spin Column	AccuPrep® Universal RNA Extraction Kit	Variable*	< 20 min
	Bacteria culture (gram-negative, gram-positive)	Spin Column	AccuPrep® Bacterial RNA Extraction Kit	Variable*	< 40 min
	Viral RNA in body fluid (blood, plasma, serum and urine)	Spin Column	AccuPrep® Viral RNA Extraction Kit	Variable*	< 30 min

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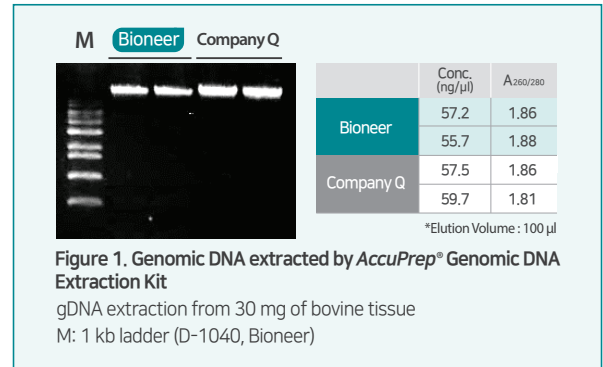
Figure 1. Genomic DNA extracted by AccuPrep® Genomic DNA Extraction Kit

gDNA extraction from 30 mg of bovine tissue
M: 1 kb ladder (D-1040, Bioneer)

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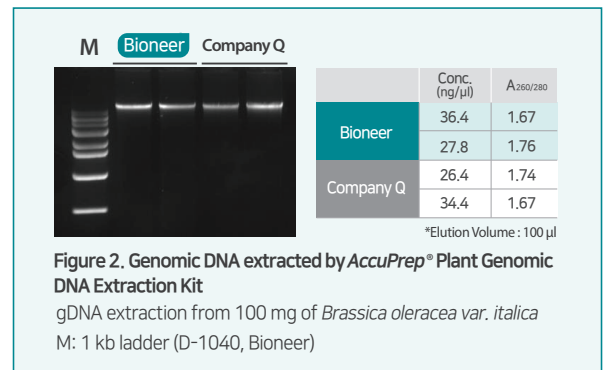
Figure 2. Genomic DNA extracted by AccuPrep® Plant Genomic DNA Extraction Kit

gDNA extraction from 100 mg of *Brassica oleracea var. italica*
M: 1 kb ladder (D-1040, Bioneer)

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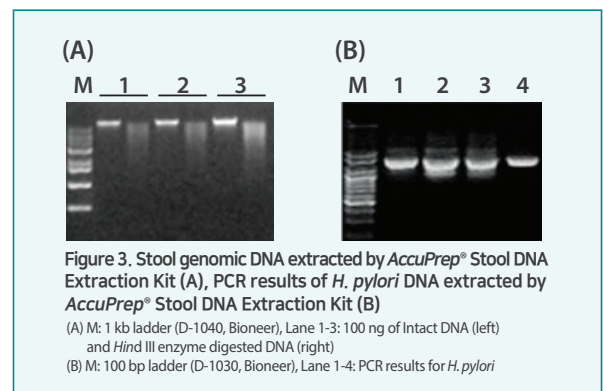
Figure 3. Stool genomic DNA extracted by AccuPrep® Stool DNA Extraction Kit (A), PCR results of *H. pylori* DNA extracted by AccuPrep® Stool DNA Extraction Kit (B)

(A) M: 1 kb ladder (D-1040, Bioneer), Lane 1-3: 100 ng of Intact DNA (left) and *Hind* III enzyme digested DNA (right)

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
Product	Genomic DNA Kit				Plant Genomic DNA Kit			Stool DNA Kit
	Whole blood	Buffy coat	Cultured cells	Mammalian tissue	Soybean	Maize	Potato	Stool
Sample	Whole blood	Buffy coat	Cultured cells	Mammalian tissue	Soybean	Maize	Potato	Stool
Expected yield	3-6 μg	15-20 μg	15-20 μg	10-15 μg	2-5 μg	1-4 μg	1-4 μg	2-5 μg
Prep. time	< 30 min				< 50 min			< 50 min
Application	Gene cloning, PCR, Real-time PCR, Southern blotting, and SNP genotyping							

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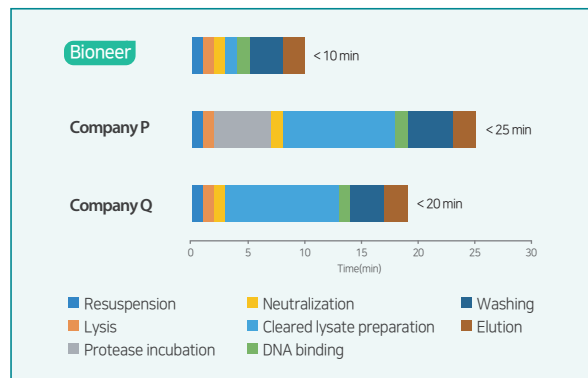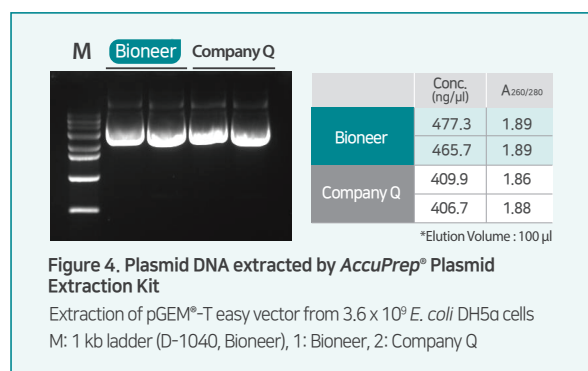
Figure 4. Plasmid DNA extracted by AccuPrep® Plasmid Extraction Kit

Extraction of pGEM®-T easy vector from 3.6×10^9 *E. coli* DH5a cells
M: 1 kb ladder (D-1040, Bioneer), 1: Bioneer, 2: Company Q

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


Product	Nano-Plus Plasmid Extraction Kit			Plasmid Mini Extraction Kit	PCR/Gel DNA Purification Kit
	Mini	Midi	Maxi		
Expected yield	Up to 30 μg	Up to 75-100 μg	Up to 300-500 μg	Up to 20 μg	Variable
Prep. time	< 10 min	< 40 min	< 60 min	< 20 min	< 30 min
Application	Sub-cloning, sequencing, transformation, transfection, and <i>in-vitro</i> transcription/translation				Sub-cloning, sequencing, labeling, and DNA concentration

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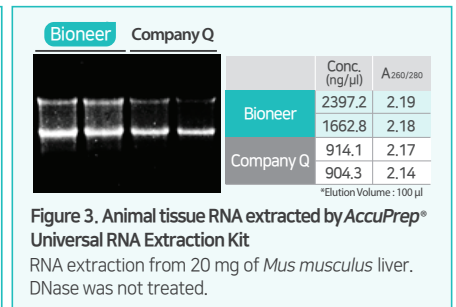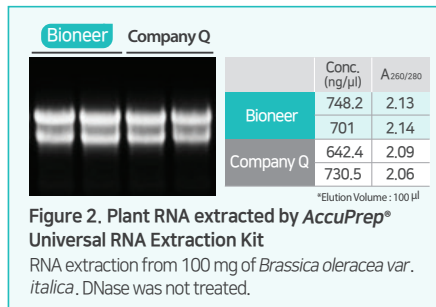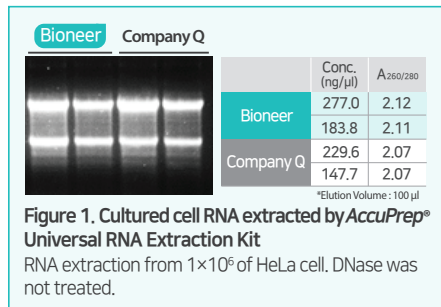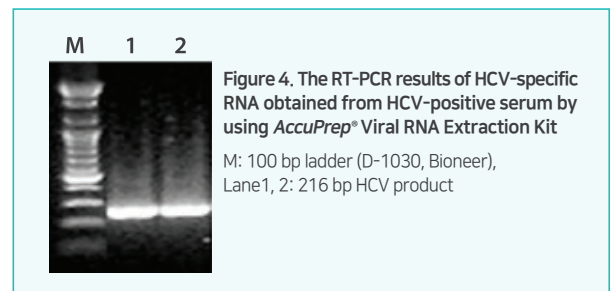
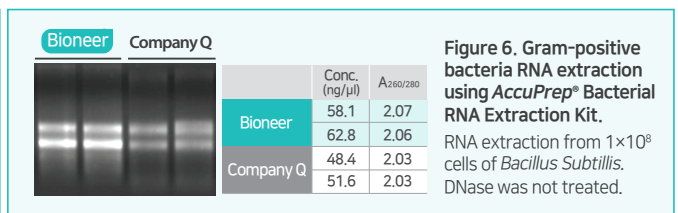
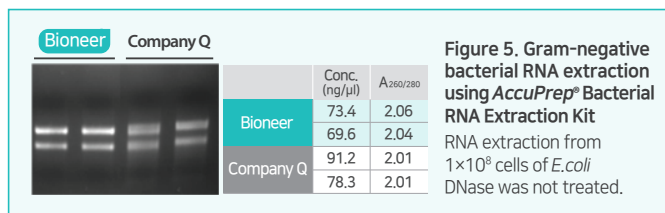
Product	Universal RNA Extraction Kit	Bacterial RNA Extraction Kit	Viral RNA Extraction Kit
Sample	Cultured animal cells, mammalian tissue, plant tissue	Bacteria culture (gram-negative, gram-positive)	Viral RNA in body fluid (blood, plasma, serum and urine)
Expected yield	Variable	Variable	Variable
Prep. time	< 20 min	< 40 min	< 30 min
Application	RT-PCR, RT-qPCR, cDNA synthesis, northern/dot/slot blot analysis, microarray, and RNA-seq		RT-PCR, RT-qPCR, cDNA synthesis, poly A+ RNA selection, northern/dot/slot blot analysis, and pathogen detection

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카탈로그 번호	제품명
K-3111/K-3112	AccuPrep® Nano-Plus Plasmid Mini Extraction Kit, 200 reactions/50 reactions
K-3122	AccuPrep® Nano-Plus Plasmid Midi Extraction Kit, 25 reactions
K-3131/K-3132	AccuPrep® Nano-Plus Plasmid Maxi Extraction Kit, 25 reactions/10 reactions
K-3030/K-3030-1	AccuPrep® Plasmid Mini Extraction Kit, 200 reactions/50 reactions
K-3032	AccuPrep® Genomic DNA Extraction Kit, 100 reactions
K-3031	AccuPrep® Plant Genomic DNA Extraction Kit, 100 reactions
K-3036	AccuPrep® Stool DNA Extraction Kit, 100 reactions
K-3140/K-3141	AccuPrep® Universal RNA Extraction Kit, 100 reactions/50 reactions
K-3142/K-3143	AccuPrep® Bacterial RNA Extraction Kit, 100 reactions/50 reactions
K-3033	AccuPrep® Viral RNA Extraction Kit, 100 reactions
K-3037/K-3038	AccuPrep® PCR/Gel DNA Purification Kit, 200 reactions/50 reactions
K-3030-2	AccuPrep® Plasmid DNA Extraction Kit for 96 well vacuum Block, 96-well plate x 2
K-3032-2	AccuPrep® Genomic DNA Extraction Kit for 96 well vacuum Block, 96-well plate x 2
K-3034-2	AccuPrep® PCR Purification Kit for 96 well vacuum Block, 96-well plate x 2

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