All about PCR

Selection Chart of Bioneer's PCR Kits

PreMix

Product	Application								
	Standard PCR	HotStart PCR	Prevent carryover contamination	PCR for gene cloning	PCR for TA cloning	High-fidelity PCR	Long range PCR	GC rich PCR	Multiplex PCR
<i>AccuPower</i> [®] PCR PreMix	√				√				
<i>AccuPower</i> [®] PCR PreMix (with UDG)	√		√		√				
<i>AccuPower</i> [®] <i>Taq</i> PCR PreMix	√				√				
<i>AccuPower</i> [®] HotStart PCR PreMix	√	√			√			√	
<i>AccuPower</i> [®] HotStart PCR PreMix (with UDG)	√	√	√		√			√	
<i>AccuPower</i> [®] <i>GoldHotStart Taq</i> PCR PreMix	√	√			√			√	
<i>AccuPower</i> [®] <i>PyroHotStart Taq</i> PCR PreMix	√	√			√			√	
<i>AccuPower</i> [®] HotStart <i>Pfu</i> PCR PreMix	√	√		√		√		√	
<i>AccuPower</i> [®] <i>ProFi Taq</i> PCR PreMix	√			√	√	√	√		
<i>AccuPower</i> [®] <i>Pfu</i> PCR PreMix	√			√		√			
<i>AccuPower</i> [®] Multiplex PCR PreMix	√	√			√				√
<i>AccuPower</i> [®] Gold Multiplex PCR PreMix	√	√			√				√

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- Conventional PCR
- Primer extension
- TA cloning
- Gene sequencing

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- Enzyme: Top DNA polymerase
- 5' → 3' exonuclease activity: No
- 3' → 5' exonuclease activity: No
- 3' - A overhang: Yes
- Fragment size: ~ 10 kb

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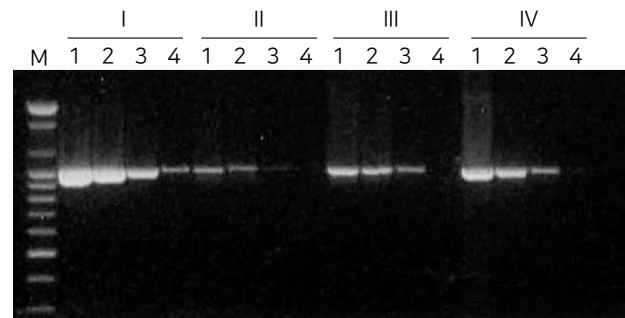
Figure 1. Comparison of sensitivity test for AccuPower® PCR PreMix and other suppliers' products using serial diluted human gDNA. Target gene: human insulin receptor gene

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- Conventional PCR
- Primer extension
- TA cloning
- Gene sequencing
- Molecular diagnosis

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- Enzyme: *Top* DNA polymerase
- 5' → 3' exonuclease activity : No
- 3' → 5' exonuclease activity : No
- 3'- A overhang : Yes
- Fragment size : ~ 10 kb

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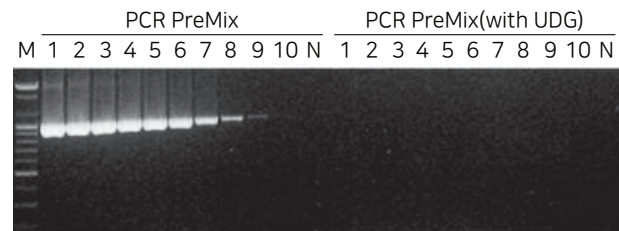
Figure 1. Efficiency of uracil DNA glycosylase using PCR products. (including uracil base)

Efficiency test of uracil DNA glycosylase was operated using serial diluted PCR products including uracil base. AccuPower® PCR PreMix was also tested for negative control. Reaction mixture was incubated at 37°C for 2 min, followed by 95°C for 5 min, 30 cycles of 20 sec at 95°C, 20 sec at 55°C, 30 sec at 72°C.

The copy number of PCR products used to test is represented below.

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- Conventional PCR
- Primer extension
- TA cloning
- Gene sequencing

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- Enzyme: Taq DNA polymerase
- 5' → 3' exonuclease activity : Yes
- 3' → 5' exonuclease activity : No
- 3'- A overhang : Yes
- Fragment size : ~ 10 kb

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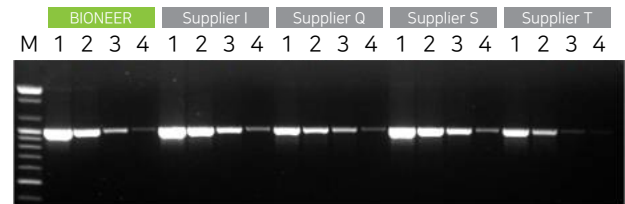
Figure 1. Comparison of PCR amplification efficiency between AccuPower® Taq PCR PreMix from Bioneer and other suppliers' PCR Master Mix.

Target gene: human insulin receptor gene.

The cycling conditions for AccuPower® Taq PCR PreMix were 95°C for 5 min, 30 cycles of 95°C for 20 sec, 55°C for 20 sec, and 72°C for 30 sec. PCR reactions using other suppliers' PCR master mix were performed according to each suppliers' protocol.

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- Primer extension
- Long range amplification from gDNA
- High amplification efficiency
- Excellent performance on difficult templates
- Amplification of low-copy targets
- High yield and high sensitivity PCR

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- Enzyme: ProFi Taq DNA polymerase
- 5' → 3' exonuclease activity: Yes
- 3' → 5' exonuclease activity: Yes
- 3' - A overhang: Yes
- Fragment size: ~ 30 kb

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Figure 1. Comparison of PCR amplification of long targets between AccuPower® ProFi Taq PCR PreMix from Bioneer and other suppliers' PCR Master Mix.

The cycling conditions for AccuPower® ProFi Taq PCR PreMix were 95°C for 5 min, 32 cycles of 95°C for 20 sec and 68°C for 15 min. PCR reactions using other suppliers' PCR master mix were performed according to each suppliers' protocol. Human gDNA was used as a template for PCR amplification.

M1: Lambda/Hind III marker (Cat. No. D-1050, Bioneer)

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- Gene synthesis
- Gene cloning
- Conventional PCR
- Primer extension
- Site-directed mutagenesis
- High fidelity PCR

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- Enzyme: Pfu DNA polymerase
- 5' → 3' exonuclease: No
- 3' → 5' exonuclease: Yes
- 3' - A overhang: No
- Fragment size: ~ 15 kb

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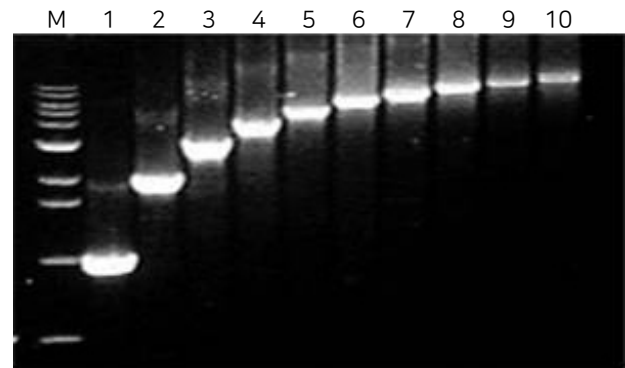
Figure 1. Amplification of lambda DNA of 1 kb to 10 kb with AccuPower® Pfu PCR PreMix.

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- Enzyme: Top DNA polymerase
- 5' → 3' exonuclease activity: No
- 3' → 5' exonuclease activity: No
- 3' - A overhang: Yes
- Fragment size: ~ 12 kb

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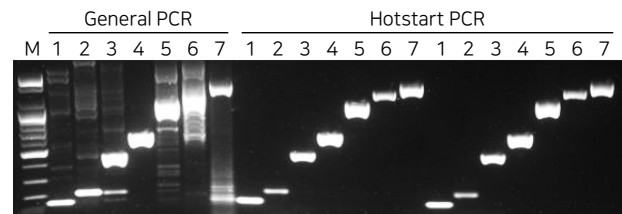
Figure 1. Specificity comparison between AccuPower® PCR PreMix and AccuPower® HotStart PCR PreMix.

- M: 100 bp DNA Ladder (Cat. No. D-1030, Bioneer)
 Lane 1: P75/P73 primer set (139 bp)
 Lane 2: P55/P53 primer set (211 bp)
 Lane 3: P55/P63 primer set (447 bp)
 Lane 4: P75/P83 primer set (618 bp)
 Lane 5: P55/P73 primer ser (1082 bp)
 Lane 6: P65/P83 primer set (1296 bp)
 Lane 7: P55/P83 primer set (1561 bp)

Figure 2. Comparison of PCR amplification sensitivity between AccuPower® HotStart PCR PreMix from Bioneer and other suppliers' HotStart PCR kit.

- M: 100 bp DNA Ladder (Cat. No. D-1030, Bioneer)
 Lane 1: Human gDNA 10 ng Lane 2: Human gDNA 1 ng
 Lane 3: Human gDNA 100 pg Lane 4: Human gDNA 10 pg

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- Enzyme: *Top* DNA polymerase
- 5' → 3' exonuclease activity: No
- 3' → 5' exonuclease activity: No
- 3' - A overhang: Yes
- Fragment size: ~ 12 kb

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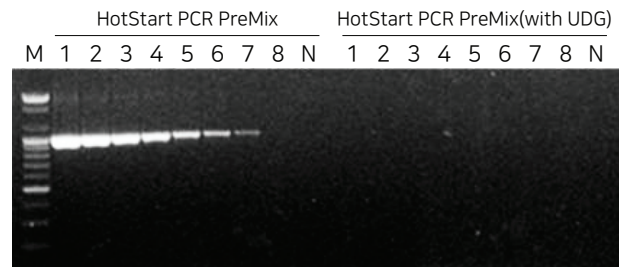
Figure 1. Efficiency of uracil DNA glycosylase using PCR product (including uracil base).

Efficiency test of uracil DNA glycosylase was operated using serial diluted PCR products including uracil base. AccuPower® HotStart PCR PreMix was also tested for negative control. Reaction mixture was incubated at 37°C for 2 min followed by 95°C for 5 min, 30 cycles of 20 sec at 95°C, 20 sec at 55°C, 30 sec at 72°C. AccuPower® HotStart PCR PreMix(with UDG) contains dUTP besides dATP, dGTP, dCTP and dTTP

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- Enzyme: Taq DNA polymerase
- 5' → 3' exonuclease activity: Yes
- 3' → 5' exonuclease activity: No
- 3' - A overhang: Yes
- Fragment size: ~ 5 kb (human)

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Figure 1. Comparison of PCR amplification specificity between AccuPower® PyroHotStart Taq PCR PreMix from Bioneer and other suppliers' HotStart PCR master mix. PCR reactions were performed according to each suppliers' protocol.

Target: Human PrP gene

The PrP gene was amplified from human gDNA with two different primer sets, separately. This data shows that AccuPower® PyroHotStart Taq PCR PreMix has higher amplification efficiency and specificity than other suppliers' HotStart PCR Master Mix.

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- Enzyme: GoldHotStart Taq DNA polymerase
- 5' → 3' exonuclease activity: Yes
- 3' → 5' exonuclease activity: No
- 3' - A overhang: Yes
- Fragment size: ~ 5 kb (human genomic DNA)

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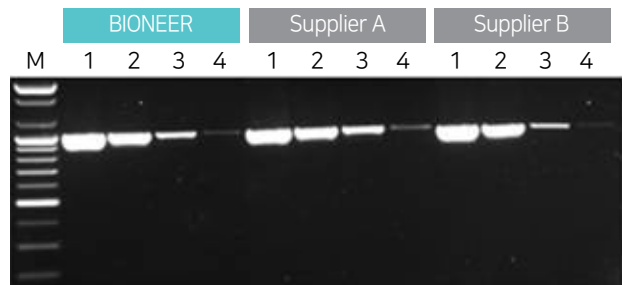
Figure 1. Comparison of PCR amplification efficiency between AccuPower® GoldHotStart Taq PCR PreMix from Bioneer and other suppliers' HotStart PCR master mix.

Target : Human insulin receptor gene.

The cycling conditions for AccuPower® GoldHotStart Taq PCR PreMix were 95°C for 5 min, 30 cycles of 95°C for 30 sec, 55°C for 30 sec and 72°C for 30 sec. PCR reactions using other suppliers' PCR master mix were performed according to each suppliers' protocol.

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- Gene cloning with blunt ends
- Site-directed mutagenesis
- High fidelity amplification
- High specificity PCR
- cDNA amplification

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- Enzyme: Pfu DNA polymerase
- 5' → 3' exonuclease activity: No
- 3' → 5' exonuclease activity: Yes
- 3' - A overhang: No
- Fragment size: ~ 5 kb

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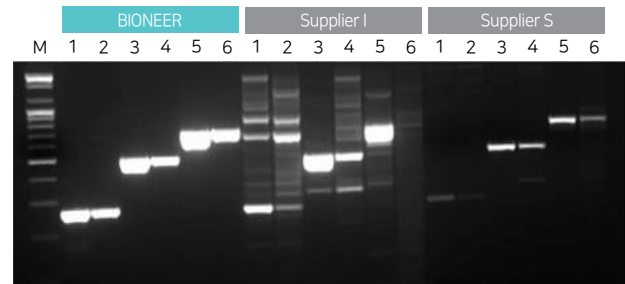
Figure 1. Comparison of PCR amplification specificity between AccuPower® HotStart Pfu PCR PreMix from Bioneer and other suppliers' Pfu DNA polymerase.

Three different primer-template systems were amplified under the same conditions. The cycling conditions were 95°C for 5 min, 35 cycles of 95°C for 30 sec, 62°C for 30 sec, and 72°C for 1 min 30 sec, and 72°C for 5 min for final extension.

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- STR analysis
- Molecular diagnostic analysis
- Genotyping assay
- Qualitative, semi-qualitative gene expression assay
- Mutant screening
- Transgenic organism analysis

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- Enzyme: HotStart Top DNA polymerase
- 5' → 3' exonuclease activity: No
- 3' → 5' exonuclease activity: No
- 3' - A overhang: Yes
- Fragment size: ~ 1 kb

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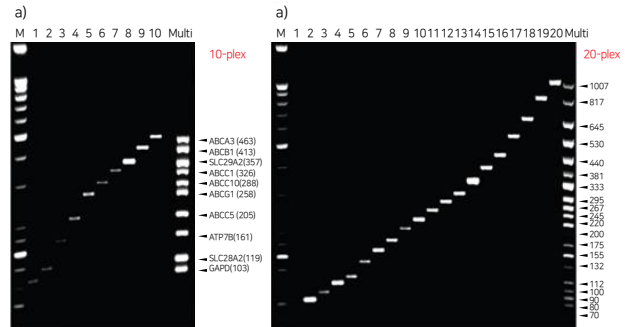
Figure 1. Single PCR and multiplex PCR using AccuPower® Multiplex PCR PreMix.

Each lane from left to right indicates the single and multiplex PCR product using AccuPower® Multiplex PCR PreMix.

a) 10-plex multiplex PCR,

b) 20-plex multiplex PCR

M: 25/100 bp Mixed DNA Ladder (Cat. No. D-1020, Bioneer)

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Target	Application
Human and Animal	STR analysis for determining genetic profiles in forensic cases
	Molecular diagnostic analysis
	Genotyping assay
	Qualitative and semi-qualitative gene expression assay
	Mutant screening
Plant	Transgenic organism analysis
	STR analysis
	Detection of pathogens/bacterial infection
	Qualitative and semi-qualitative gene expression assay

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- Enzyme: Top DNA polymerase
- 5' → 3' exonuclease activity: No
- 3' → 5' exonuclease activity: No
- 3' - A overhang: Yes
- Fragment size: ~ 1 kb

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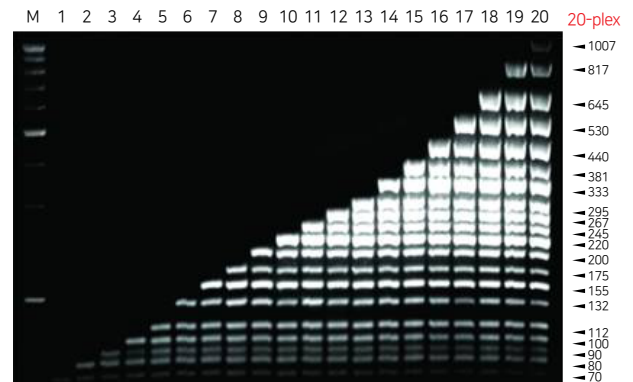
Figure 1. High specificity of AccuPower® Gold Multiplex PCR PreMix. Each lane from left to right represents the progressive number of primer sets (1 - 20) included in single AccuPower® Gold Multiplex PCR PreMix reactions.

Rxn. Condition : 95°C for 10 min, followed by 35 cycles of 30 sec at 95°C, 30 sec at 57°C, 60 sec at 72°C

M: 25/100 bp Mixed DNA Ladder (Cat. No. D-1020, Bioneer)

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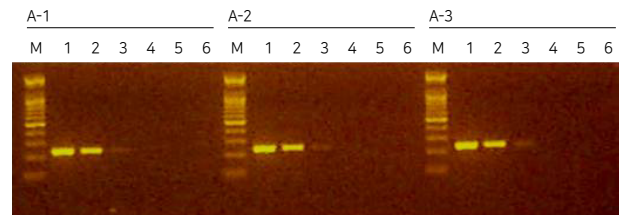
Figure 1. Custom AccuPower® PCR PreMix: Reproducibility & Sensitivity. Custom AccuPower® DsRed2 PCR PreMix was tested three times, A-1, A-2 and A-3, with the condition of: 1. 94°C 2 min; 2. 30 cycles of 94°C 30 sec, 58°C 30 sec and 72°C 30 sec; 3. 72°C 3 min. 5 ul of amplified PCR product was run on a 2% TBE agarose gel.

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PreMix Type

카탈로그 번호	제품명
<i>AccuPower</i> [®] PCR PreMix	
K-2012 / K-2013	<i>AccuPower</i> [®] PCR PreMix, 96 tubes, 20 µl / 50 µl
K-2016 / K-2017	<i>AccuPower</i> [®] PCR PreMix, 480 tubes, 20 µl / 50 µl
K-2011	<i>AccuPower</i> [®] PCR PreMix, 0.5 ml thin-wall tube with attached cap, 100 tubes, 50 µl
<i>AccuPower</i> [®] PCR PreMix (with UDG)	
K-2012-1	<i>AccuPower</i> [®] PCR PreMix (with UDG), 96 tubes, 20 µl
K-2016-1	<i>AccuPower</i> [®] PCR PreMix (with UDG), 480 tubes, 20 µl
<i>AccuPower</i> [®] Taq PCR PreMix	
K-2601 / K-2603	<i>AccuPower</i> [®] Taq PCR PreMix, 96 tubes, 20 µl / 50 µl
K-2602 / K-2604	<i>AccuPower</i> [®] Taq PCR PreMix, 480 tubes, 20 µl / 50 µl
<i>AccuPower</i> [®] ProFi Taq PCR PreMix	
K-2631 / K-2633	<i>AccuPower</i> [®] ProFi Taq PCR PreMix, 96 tubes, 20 µl / 50 µl
K-2632 / K-2634	<i>AccuPower</i> [®] ProFi Taq PCR PreMix, 480 tubes, 20 µl / 50 µl
<i>AccuPower</i> [®] Pfu PCR PreMix	
K-2022 / K-2023	<i>AccuPower</i> [®] Pfu PCR PreMix, 96 tubes, 20 µl / 50 µl
K-2024 / K-2025	<i>AccuPower</i> [®] Pfu PCR PreMix, 480 tubes, 20 µl / 50 µl
K-2027	<i>AccuPower</i> [®] Pfu PCR PreMix, 0.5 ml thin-wall tubes with attached cap, 100 tubes, 50 µl
<i>AccuPower</i> [®] HotStart PCR PreMix	
K-5050 / K-5052	<i>AccuPower</i> [®] HotStart PCR PreMix, 96 tubes, 20 µl / 50 µl
K-5051 / K-5057	<i>AccuPower</i> [®] HotStart PCR PreMix, 480 tubes, 20 µl / 50 µl
K-5053	<i>AccuPower</i> [®] HotStart PCR PreMix, 0.5 ml thin-wall tubes with attached cap, 100 tubes, 50 µl
<i>AccuPower</i> [®] HotStart PCR PreMix (with UDG)	
K-5050-1	<i>AccuPower</i> [®] HotStart PCR PreMix (with UDG), 96 tubes, 20 µl
K-5051-1	<i>AccuPower</i> [®] HotStart PCR PreMix (with UDG), 480 tubes, 20 µl
<i>AccuPower</i> [®] PyroHotStart Taq PCR PreMix	
K-2611 / K-2613	<i>AccuPower</i> [®] PyroHotStart Taq PCR PreMix, 96 tubes, 20 µl / 50 µl
K-2612 / K-2614	<i>AccuPower</i> [®] PyroHotStart Taq PCR PreMix, 480 tubes, 20 µl / 50 µl
<i>AccuPower</i> [®] GoldHotStart Taq PCR PreMix	
K-2621 / K-2623	<i>AccuPower</i> [®] GoldHotStart Taq PCR PreMix, 96 tubes, 20 µl / 50 µl
K-2622 / K-2624	<i>AccuPower</i> [®] GoldHotStart Taq PCR PreMix, 480 tubes, 20 µl / 50 µl
<i>AccuPower</i> [®] HotStart Pfu PCR PreMix	
K-2301 / K-2302	<i>AccuPower</i> [®] HotStart Pfu PCR PreMix, 96 tubes, 20 µl / 50 µl
K-2303 / K-2304	<i>AccuPower</i> [®] HotStart Pfu PCR PreMix, 480 tubes, 20 µl / 50 µl
<i>AccuPower</i> [®] Multiplex PCR PreMix	
K-2111 / K-2112	<i>AccuPower</i> [®] Multiplex PCR PreMix, 96 tubes, 20 µl / 50 µl
K-2113 / K-2114	<i>AccuPower</i> [®] Multiplex PCR PreMix, 480 tubes, 20 µl / 50 µl
<i>AccuPower</i> [®] Gold Multiplex PCR PreMix	
K-2115 / K-2117	<i>AccuPower</i> [®] Gold Multiplex PCR PreMix, 96 tubes, 20 µl / 50 µl
K-2116 / K-2118	<i>AccuPower</i> [®] Gold Multiplex PCR PreMix, 480 tubes, 20 µl / 50 µl

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